

Agilent Cary UV-Vis (NIR) Spectrophotometers Models 4000, 5000, 6000i and 7000

Long Term Shut Down and Start Up Instructions

Shut down the system

To shut down the system:

- 1 Turn off the instrument purge from the supply (if applicable).
- 2 Turn off the instrument power from the panel switch.
- 3 Disconnect the mains power lead.
- 4 Remove any installed accessories and store them in their original packaging.
- 5 Fit the sample compartment door.
- 6 Close the sample compartment lid.
- 7 Shutdown the controlling PC.

Start up the system

To start up the system:

- 1 Turn on the instrument purge from the supply (if applicable).
- 2 Connect the instrument power lead.
- 3 Switch on the instrument.
- 4 Switch on the controlling PC.
- 5 Allow a 2-hour instrument warmup period.
- 6 Open the Validate application and run the Instrument Performance Tests.