

Analyzing Chemical Composition of Rigid Polyurethane (PU) Foams using FTIR Microscopy and DRIFTS-IR

Pittcon 2015 2250-4P

John Kinyanjui, Ph.D., Jeff Head, Mark Talbott, Ph.D.;
Shimadzu Scientific Instruments, Columbia, MD, USA

Analyzing Chemical Composition of Rigid Polyurethane (PU) Foams using FTIR Microscopy and DRIFTS-IR

Introduction


- The chemical composition of rigid polyurethane foams was studied using FTIR microscopy and diffuse reflectance Fourier transform spectroscopy (DRIFTS) FTIR.
- Temperature gradients during the processing of rigid polyurethane (PU) foams leads to chemical gradients that might be observed across the radial and axial dimensions of the foam.
- Unreacted reactants and products from side reactions can remain trapped in the cavities of the foam hampering mechanical strength.
- A highly **complex and heterogeneous** polymeric composite results.
- The chemical composition of the cellular structure of the foam was studied with distinct spectra of the gas filled cell nucleus and cell wall.
- FTIR spectroscopy offers an ideal tool for studying variations in chemical composition at both bulk and microscopic levels.

Experimental Conditions - Reactants


- Commercially available rigid polyurethane pour foam starting materials (designated as Part A and PART B by manufacturer) were used.
- Although the exact formulation was not provided
 - PART A normally contains the diisocyanate.
 - Part B normally contains the polyol, water (blowing agent), a catalyst and a surfactant premixed at optimized ratios.
- Equal amounts of PART A and PART B were mixed and stirred until the mixture appeared uniform.
- The foam was then allowed to rise and "cure" over a period of five (5) days.
- Foam was sectioned and sample measurement carried out on the fifth day.
- Based on FTIR results unique **uretoneimine** crosslinking functionality might be present in the foam giving clues on exact formulation of starting materials. ^[1,2]

Rigid PU Foam Formation Reaction

General Reaction ^[1,2]

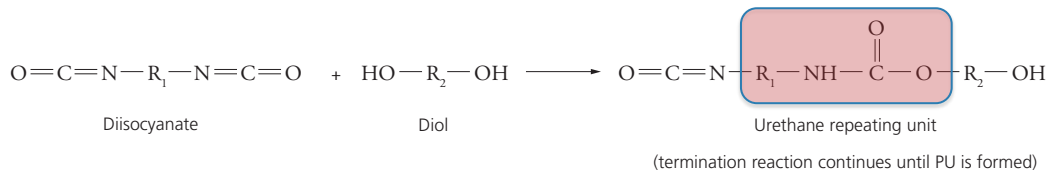


Foam Blowing Reaction ^[1,2]

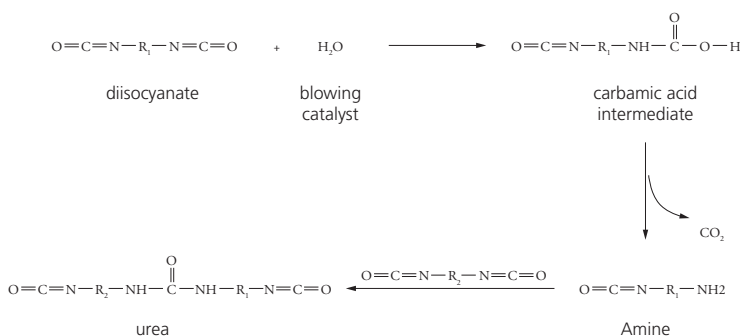


Analyzing Chemical Composition of Rigid Polyurethane (PU) Foams using FTIR Microscopy and DRIFTS-IR

Experimental Conditions - Foam Mold Designation

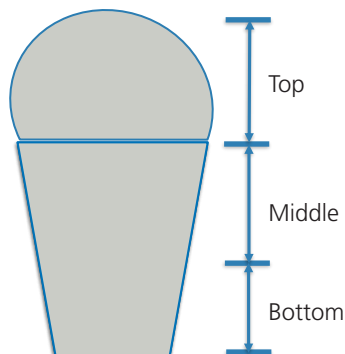

- Three sections in the mold were cut and designated as "top", "middle" and "bottom" for sampling.
- Samples were bored out from the radial center of each area.
- For the middle section, three samples were bored out and designated as "center", "Edge 1" and "Edge 2".

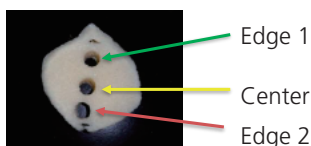

Instrumentation



The AIM-8800 connected to IRTracer-100



CO₂ filled cell cell wall
Light Microscope Image of PU Foam

Sample Preparation

- Sample bored out from the center of the middle section of the mold was tested. A cylinder microtome was used to produce thin specimens for microscopy.




The PIKE EasiDiff DRIFTS Accessory

Sample Preparation

- Pristine samples were bored out and sliced to fit tightly into the 10 mm diameter sample cup provided with the accessory.


Analyzing Chemical Composition of Rigid Polyurethane (PU) Foams using FTIR Microscopy and DRIFTS-IR

DRIFTS-IR: Top to Bottom of Mold

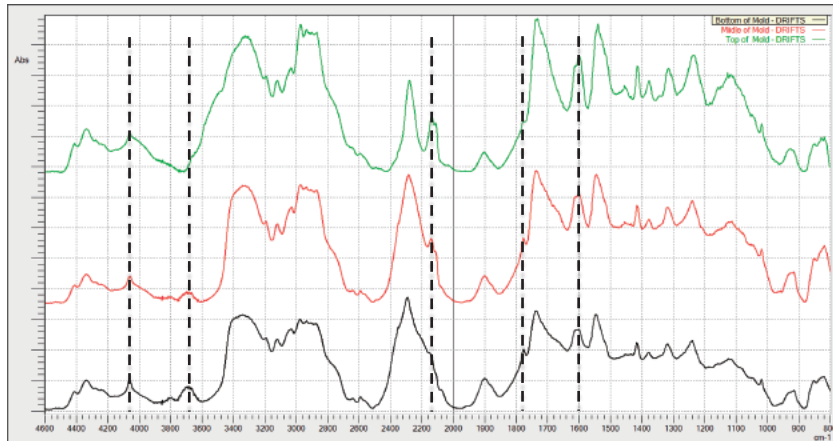


Composition of the mold changes as you move from top to bottom of the mold.

DRIFTS IR: Radial Profile of Middle Section of Mold

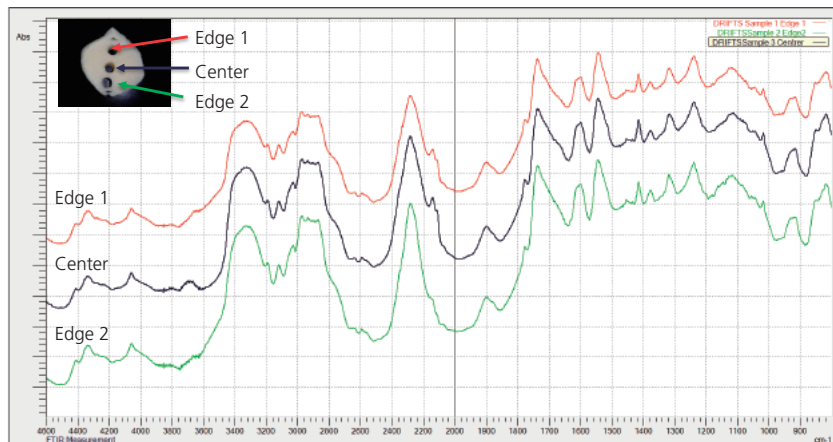

Composition across the radial axis of the middle section does not change appreciably.


FTIR Microscopy: Cell and Cell Wall Spectra

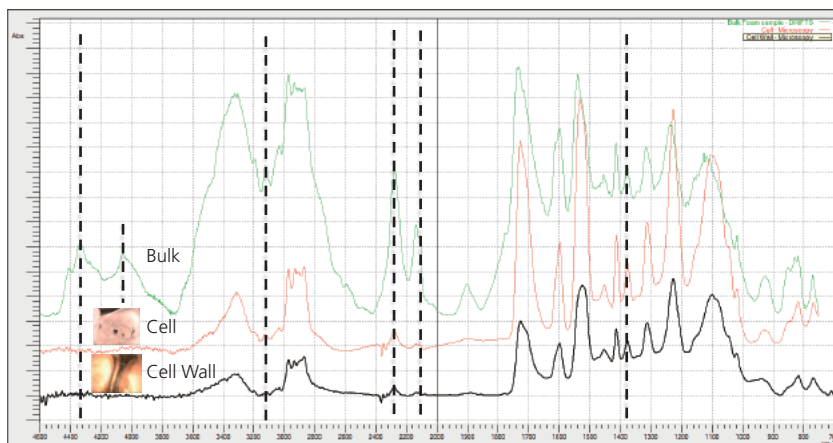

Composition of cell and cell wall different from the bulk foam due to trapped reactants and side reaction products (such as urea) in foam micro-pores.

Analyzing Chemical Composition of Rigid Polyurethane (PU) Foams using FTIR Microscopy and DRIFTS-IR


Cross-Linking in PU Foam^[1,2]

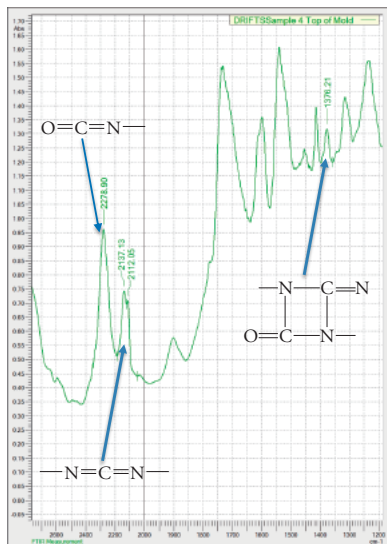


Unreacted Isocyanate reacts by itself to form *carbodiimide*

Carbodiimide reacts with isocyanate to form *urethaneimine* crosslinks

Urethaneimine linkages lead to increased rigidity

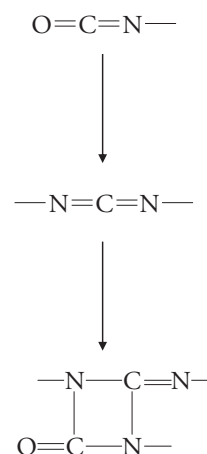

DRIFTS-IR: Top to Bottom of Mold

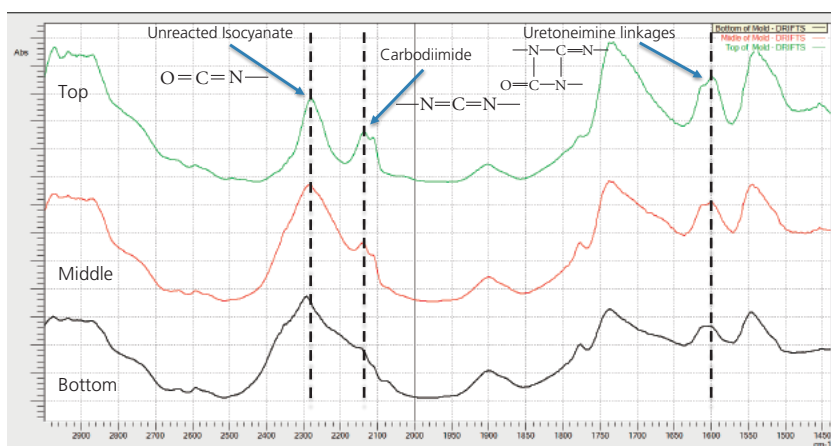

Composition of the mold changes as you move from top to bottom of the mold.

Conclusions

- FTIR microscopy was successfully used to show distinct chemical composition at the cellular level as compared to the bulk composition.
- With DRIFTS-IR the spectra of pristine foam samples can be easily collected with minimal sample preparation to provide the bulk chemical properties of the foam.
- Foam aging and curing including the level of cross-linking can also be monitored using DRIFTS-IR.
- The mechanical properties of the PU foam is highly dependent on the chemical properties of the foam.^[1,2] FTIR data can be used as a tool to assist in tuning the mechanical properties of the foam.
- Previous work has shown that the level of cross-linking in PU foam can be directly correlated to the Young's modulus of the foam.^[1,2]

Analyzing Chemical Composition of Rigid Polyurethane (PU) Foams using FTIR Microscopy and DRIFTS-IR

References

1. Mohan, R.B.; O'Toole B.J.; Malpica, J.; Hatchett D.W.; Kodippili G.; and Kinyanjui J.M. **Effects of Processing Temperature on ReCrete Polyurethane Foam.** Journal of Cellular Plastics 2008; 44: 327-345.
2. Hatchett, D.W.; Kodippili, G; Kinyanjui, J.M.; Benincasa, F.; Sapochak, L. **FTIR Analysis of Thermally Processed PU Foam.** Polymer Degradation and Stability; 2005; 87(3), 555-561.